# Set Your Sights on Superior Performance

## PAL Smart SPME Arrow

- Rugged, stainless-steel construction ensures longer lifetimes.
- Faster extraction means higher sample throughput.
- Better sensitivity allows lower LODs.
- Now with Smart technology.



Pure Chromatography



## **Set Your Sights on Superior Performance**

Solid phase microextraction (SPME) is a fast, automated sample preparation technology that reduces sample handling, extraction time, and solvent consumption, making it a popular choice in environmental, food, and clinical laboratories. However, traditional SPME fiber technology has some significant drawbacks, including poor mechanical stability and a small phase volume.

The PAL Smart SPME Arrow system is a revolutionary change in microextraction that combines exceptional robustness with faster extraction times and trace-level sensitivity (Table I). In contrast to traditional SPME fibers, SPME Arrows contain significantly more phase volume, which allows more target analyte to be extracted in less time. In addition, the stainless-steel construction, unique Arrow tip, inner stabilizing rod, and outer sheath design fully protect the phase, minimizing both mechanical damage and analyte loss during sample transfer.



Table I: Comparing Restek PAL SPME Arrow to Traditional SPME Fiber (Headspace Technique)

	Traditional Fibers	1.1 mm Arrow	1.5 mm Arrow	The Arrow Advantage
Rugged, stainless-steel construction	No	Yes	Yes	<ul> <li>Longer lifetimes: Arrow construction protects the phase and minimizes both damage and analyte loss.</li> </ul>
• Protective sheath				• Less downtime due to breakage.
<ul> <li>Stabilizing inner rod</li> </ul>				• Eliminates septa coring.
Arrow-shaped tip				
Surface area	9.4 mm <sup>2</sup>	44 mm <sup>2</sup>	63 mm <sup>2</sup>	Increase sample throughput—higher surface area lets you reduce extraction time and <b>analyze more samples per shift.</b>
Phase (PDMS) volume	0.6 µL	3.8 µL	12 µL	With significantly more phase, SPME Arrow provides better sensitivity so you can <b>lower detection limits with confidence.</b>

Our product line is continually expanding! See what's new at www.restek.com/SPME



## **On Target: Longer Lifetime**

One of the main drawbacks to SPME fibers is their fragility. Even during routine use, they are easily damaged and can be irreparably broken (Figure 1). Busy labs will benefit greatly from the reliability of rugged Restek PAL Smart SPME Arrow. The unique, stainless-steel Arrow design includes a robust inner stabilizing rod that resists bending and breakage as well as an outer sheath that protects the phase coating and prevents both physical damage and analyte loss. Typically, SPME Arrows perform well for hundreds of extractions, lasting two to three times longer than traditional SPME fibers. In addition, the unique arrow-shaped tip pierces the septum cleanly and with little resistance, extending septum lifetime (Figure 2).

#### Figure 1: SPME fibers break easily during routine operation.



**Figure 2:** The rugged construction of the Restek PAL SPME Arrow prevents breakage under normal use.



### Which Restek PAL SPME Arrow is best for my application?

Restek PAL Smart SPME Arrows are suitable for a wide range of analyte chemistries and sample matrices. They are suitable for manual injection and are fully backwards compatible with any generation of PAL3 Systems. Choose the best SPME Arrow for your application based on the properties of your target compounds.

- Trace analysis in foodstuffs
- Drugs and pharmaceuticals
- Herbicides/pesticides

- Medical diagnostics
- Trace impurities in polymers and solid samples
- · Solvent residues in raw materials
- Water analysis (organics in water)

Check your instrument manufacturer requirements or contact Restek Technical Support at **www.restek.com/contact-us** for further assistance.



## **On Target: Higher Sample Throughput**

Restek PAL Smart SPME Arrows provide much faster extraction times than traditional SPME fibers because Arrows have much more phase volume. As shown in Figure 3, the increased phase volume allows more target analyte to be extracted in just seconds using an Arrow than can be extracted from a traditional fiber. Being able to extract what you need in a fraction of the time means more samples can be analyzed per day, which improves lab efficiency and profitability. The example in Table II demonstrates a nearly 50% increase in productivity!

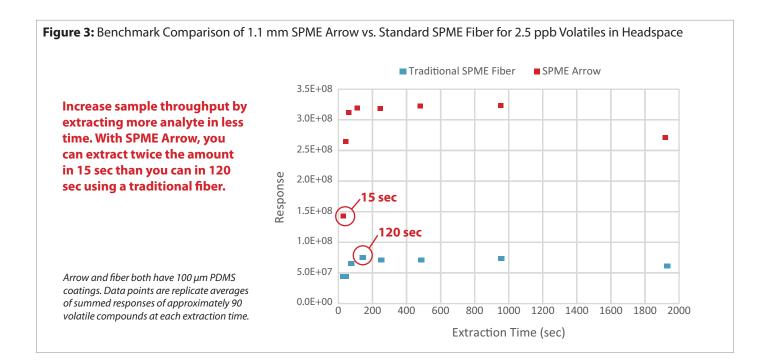


 Table II: Analyze more samples per day with Restek PAL SPME Arrow.

	Sample Collection and Desorption Steps with Associated Times (sec)								
Target Analytes	Select Sample Vial	Vial Incubation / Fiber Conditioning	Vial Transfer	Sample Extraction (Headspace)	GC Equilibration	Desorb Fiber	Return Sample Vial	Total Time	Samples per Day
Traditional SPME Fiber	22	120	20	120	15	10	15	322	268
SPME Arrow	22	120	20	15	15	10	15	217	<b>398</b> (~50% increase)

Note: Processes that take <10 sec to perform were omitted from the table. Actual samples per day is dependent on GC cycle time.






## **On Target: Lower Detection Limits**

Developing a new method or trying to improve the performance of an existing one? Restek PAL Smart SPME Arrows have more phase volume, so they provide much greater sensitivity than traditional fibers. To demonstrate this, 82 volatile compounds were analyzed under three different headspace extraction time and extraction volume combinations. Results under all experimental conditions definitively show that Arrow outperforms traditional SPME fiber and provides much higher analyte responses (Table III and Figure 4).


Table III: Analyte response is much higher with SPME Arrow than with a traditional SPME fiber.

		Average % Increase in	n Response of Arrow vs. Traditional Fiber
Extraction Time (min)	Extraction Volume (mL Water)	1.1 mm Arrow	1.5 mm Arrow
10	10	297%	527%
5	10	618%	896%
10	5	446%	634%

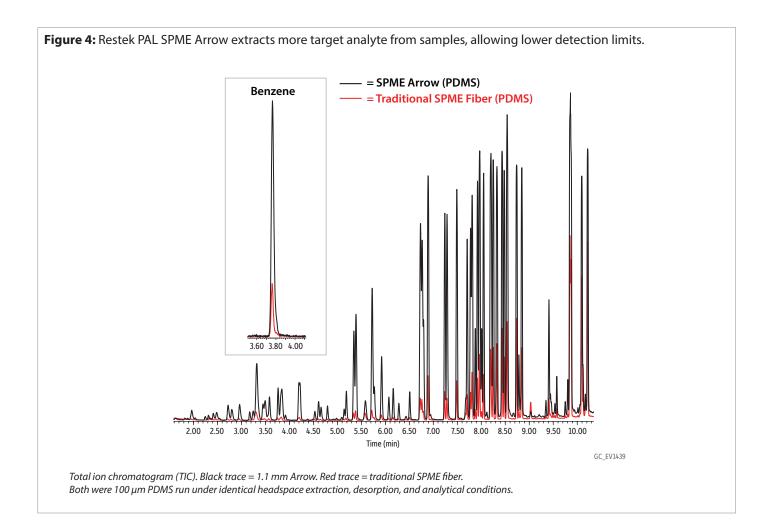




Turning to the emerging cannabis market and the triple-phase (DVB/Carbon WR/PDMS) Arrow, you can also expect significantly higher analyte responses with residual solvents and other classes of compound (Figure 5).

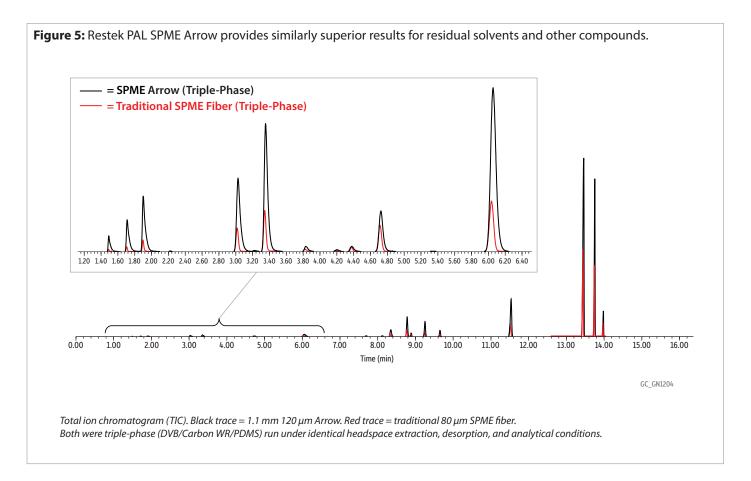

Even after 70 minutes of immersion extraction, the compound concentration of PAHs for the traditional SPME fiber could not equal the amount obtained using a SPME Arrow (Figure 6).

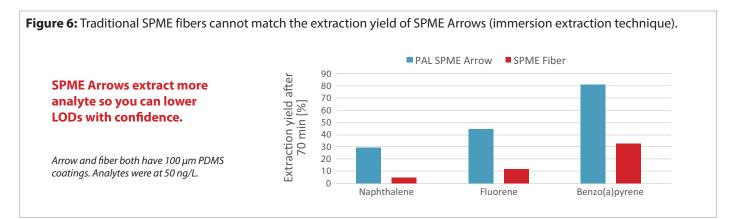

Switch to SPME Smart Arrow, and Set Your Sights on Superior Performance!



## **Get set up for Superior SPME!**

#### **Restek PAL Smart SPME Arrow**

- Rugged, stainless-steel construction ensures longer lifetimes.
- Faster extraction means higher sample throughput.
- Better sensitivity allows lower LODs.
- Choose format based on sampling technique:
  - •1.1 mm sleeve, if only using headspace sampling.
  - 1.5 mm wide sleeve, if using direct immersion sampling (works for headspace, too).
- Now with Smart technology—each SPME Arrow is equipped with a unique Smart chip containing parameters, ranges, and usage history.

Recommended maximum GC inlet pressure is 50 psi or less.

All PAL SPME Smart Arrows have 20 mm of phase bonded onto stainless steel.

**PAL system compatibility:** Smart SPME Arrows are fully backwards compatible with any generation of PAL3 Systems. Additional hardware will be also required, including the SPME Arrow Tool, Heatex Stirrer, Arrow Conditioning Module, and Agitator Module with the perforated lid. The SPME Arrow is not supported on the HTS/Combi PAL systems.



Due to the relatively large diameter of PAL Smart SPME Arrows, you must modify the GC inlet using an instrument-specific conversion kit from Restek prior to use. Converted inlets are compatible with all standard injection techniques (SPME Arrow, liquid syringe, headspace, etc.); no need to switch inlets after installation.

Product Name	Color	Diameter	Material	<b>Recommended Analytes</b>	Units
PAL Smart SPME Arrow Method Development Kit					kit
PAL Smart SPME Arrow 1.10 mm: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, ea.	Light Blue	1.1 mm	120 µm Carbon-WR/PDMS (Carbon Wide Range/ Polydimethylsiloxane)	Highly volatile, 30–225 g/mol*	ea.
PAL Smart SPME Arrow 1.10 mm: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, 3-pk.	Light Blue	1.1 mm	120 µm Carbon-WR/PDMS (Carbon Wide Range/ Polydimethylsiloxane)	Highly volatile, 30–225 g/mol*	3-pk.
PAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, ea.	Dark Gray	1.1 mm	120 µm DVB/Carbon-WR/PDMS (Divinylbenzene/ Carbon Wide Range/Polydimethylsiloxane)	Volatile and semivolatile, 40–275 g/mol*	ea.
PAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, 3-pk.	Dark Gray	1.1 mm	120 µm DVB/Carbon-WR/PDMS (Divinylbenzene/ Carbon Wide Range/Polydimethylsiloxane)	Volatile and semivolatile, 40–275 g/mol*	3-pk.
PAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, ea.	Violet	1.1 mm	120 µm DVB/PDMS (Divinylbenzene/ Polydimethylsiloxane)	Amines and polar compounds, 60–300 g/mol*	ea.
PAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, 3-pk.	Violet	1.1 mm	120 µm DVB/PDMS (Divinylbenzene/ Polydimethylsiloxane)	Amines and polar compounds, 60–300 g/mol*	3-pk.
PAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, ea.	Red	1.1 mm	100 µm PDMS (Polydimethylsiloxane)	Volatile, 60–275 g/mol*	ea.
PAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, 3-pk.	Red	1.1 mm	100 µm PDMS (Polydimethylsiloxane)	Volatile, 60–275 g/mol*	3-pk.
PAL Smart SPME Arrow 1.10 mm: PA, Phase Thickness 100 um, Phase Length 20 mm, Gray, ea.	Gray	1.1 mm	100 µm PA (Polyacrylate)	Polar, semivolatile, 80–300 g/mol*	ea.
PAL Smart SPME Arrow 1.10 mm: PA, Phase Thickness 100 um, Phase Length 20 mm, Gray, 3-pk	Gray	1.1 mm	100 µm PA (Polyacrylate)	Polar, semivolatile, 80–300 g/mol*	3-pk.
PAL Smart SPME Arrow 1.5 mm: PDMS, Phase Thickness 250 um, Phase Length 20 mm, Black, ea.	Black	1.5 mm	250 µm PDMS (Polydimethylsiloxane)	Volatile, 60–275 g/mol (high capacity)*	ea.
PAL Smart SPME Arrow 1.5 mm: PDMS, Phase Thickness 250 um, Phase Length 20 mm, Black, 3-pk.	Black	1.5 mm	250 µm PDMS (Polydimethylsiloxane)	Volatile, 60–275 g/mol (high capacity)*	3-pk.
PAL Smart SPME Arrow 1.50 mm Wide Sleeve: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, ea.	Light Blue	1.5 mm wide sleeve	120 µm Carbon-WR/PDMS (Carbon Wide Range/ Polydimethylsiloxane)	Highly volatile, 30–225 g/mol*	ea.
PAL Smart SPME Arrow 1.50 mm Wide Sleeve: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, 3-pk.	Light Blue	1.5 mm wide sleeve	120 µm Carbon-WR/PDMS (Carbon Wide Range/ Polydimethylsiloxane)	Highly volatile, 30–225 g/mol*	3-pk.
PAL Smart SPME Arrow 1.50 mm Wide Sleeve: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, ea.	Dark Gray	1.5 mm wide sleeve	120 µm DVB/Carbon-WR/PDMS (Divinylbenzene/ Carbon Wide Range/Polydimethylsiloxane)	Volatile and semivolatile, 40–275 g/mol*	ea.
PAL Smart SPME Arrow 1.50 mm Wide Sleeve: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, 3-pk.	Dark Gray	1.5 mm wide sleeve	120 µm DVB/Carbon-WR/PDMS (Divinylbenzene/ Carbon Wide Range/Polydimethylsiloxane)	Volatile and semivolatile, 40–275 g/mol*	3-pk.
PAL Smart SPME Arrow 1.50 mm Wide Sleeve: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, ea.	Violet	1.5 mm wide sleeve	120 µm DVB/PDMS (Divinylbenzene/ Polydimethylsiloxane)	Aromatic semivolatile, 60–300 g/mol*	ea.
PAL Smart SPME Arrow 1.50 mm Wide Sleeve: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, 3-pk.	Violet	1.5 mm wide sleeve	120 µm DVB/PDMS (Divinylbenzene/ Polydimethylsiloxane)	Aromatic semivolatile, 60–300 g/mol*	3-pk.
PAL Smart SPME Arrow 1.50 mm Wide Sleeve: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, ea.	Red	1.5 mm wide sleeve	100 µm PDMS (Polydimethylsiloxane)	Volatile, 60–275 g/mol*	ea.
PAL Smart SPME Arrow 1.50 mm Wide Sleeve: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, 3-pk.	Red	1.5 mm wide sleeve	100 µm PDMS (Polydimethylsiloxane)	Volatile, 60–275 g/mol*	3-pk.
	<ul> <li>PAL Smart SPME Arrow Method Development Kit</li> <li>PAL Smart SPME Arrow 1.10 mm: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, ea.</li> <li>PAL Smart SPME Arrow 1.10 mm: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, ea.</li> <li>PAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, ea.</li> <li>PAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, 3-pk.</li> <li>PAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, ea.</li> <li>PAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, a-pk.</li> <li>PAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Violet, 3-pk.</li> <li>PAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, ea.</li> <li>PAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, a-pk.</li> <li>PAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Gray, ea.</li> <li>PAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Gray, a-pk.</li> <li>PAL Smart SPME Arrow 1.10 mm: PA, Phase Thickness 100 um, Phase Length 20 mm, Gray, a-pk.</li> <li>PAL Smart SPME Arrow 1.50 mm: PDMS, Phase Thickness 250 um, Phase Length 20 mm, Black, a-pk.</li> <li>PAL Smart SPME Arrow 1.50 mm Wide Sleeve: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, a-Ph.</li> <li>PAL Smart SPME Arrow 1.50 mm Wide Sleeve: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, 3-pk.</li> <li>PAL Smart SPME Arrow 1.50 mm Wide Sleeve: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, a-pk.</li> <li>PAL Smart SPME Arrow 1.50 mm Wide Sleeve: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, a-pk.</li> <li>PAL Smart SPME Arrow 1.50 mm Wide Sleeve: DVB/Carbon-WR/PDMS, Phase Thickness 120</li></ul>	PAL Smart SPME Arrow Method Development KitLight BluePAL Smart SPME Arrow 1.10 mm: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, a.e.Light BluePAL Smart SPME Arrow 1.10 mm: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, 3-pk.Dark GrayPAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, ea.Dark GrayPAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Dark Gray, 3-pk.Dark GrayPAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, ea.VioletPAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, 3-pk.VioletPAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, a-pk.RedPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, 3-pk.RedPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Gray, a-pk.GrayPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Gray, a-pk.GrayPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 250 um, Phase Length 20 mm, Gray, a-pk.BlackPAL Smart SPME Arrow 1.5 mm: PDMS, Phase Thickness 250 um, Phase Length 20 mm, Black, ea.Light BluePAL Smart SPME Arrow 1.5 mm: PDMS, Phase Thickness 250 um, Phase Length 20 mm, Black, a-pk.Light BluePAL Smart SPME Arrow 1.50 mm Wide Sleeve: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, a-pk.Dark GrayPAL Smart SPME Arrow 1.50 mm Wide Sleeve	PAL Smart SPME Arrow Method Development KitLight Blue1.1 mmPAL Smart SPME Arrow 1.10 mm: Carbon-WR/PDMS, Phase Thickness 120 um, PAL Smart SPME Arrow 1.10 mm: Carbon-WR/PDMS, Phase ThicknessLight Blue1.1 mmPAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase ThicknessDark Gray1.1 mmPAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase ThicknessDark Gray1.1 mmPAL Smart SPME Arrow 1.10 mm: DVB/Carbon-WR/PDMS, Phase ThicknessDark Gray1.1 mmPAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase ThicknessDark Gray1.1 mmPAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 120 um, PhaseLength 20 mm, Violet, ea.Violet1.1 mmPAL Smart SPME Arrow 1.10 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, a.Red1.1 mmPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, a.Red1.1 mmPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, a.Red1.1 mmPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, a.Gray1.1 mmPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Gray, ea.Gray1.1 mmPAL Smart SPME Arrow 1.10 mm: PDMS, Phase Thickness 200 um, Phase Length 20 mm, Red, a.SimmSimmPAL Smart SPME Arrow 1.50 mm Wide Sleeve: Carbon-WR/PDMS, PhaseBlack1.5 mmPAL Smart SPME Arrow 1.50 mm Wide Sleeve: Carbon-WR/PDMS, PhaseLight BlueSimmPAL Smart SPME Arrow 1.50 mm Wide Sleeve: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Le	PAL Smart SPME Arrow Method Development Kit       120 µm Carbon-WR/PDMS (Carbon Wide Range/ Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: Carbon-WR/PDMS, Phase Thickness 120 µm, PAL Smart SPME Arrow 110 mm: Carbon-WR/PDMS, Phase Thickness       11 µm       120 µm Carbon-WR/PDMS (Carbon Wide Range/ Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: Carbon-WR/PDMS, Phase Thickness       Dark Gray       1.1 µm       120 µm DVB(Carbon-WR/PDMS, Phase Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: DVB(Carbon-WR/PDMS, Phase Thickness       Dark Gray       1.1 µm       120 µm DVB(Carbon-WR/PDMS, Phase Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: DVB(PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, a.       1.1 µm       120 µm DVB(PDMS (Divinylbenzene/ Carbon Wide Range/Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Violet, a.       1.1 µm       120 µm DVB/PDMS (Divinylbenzene/ Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, ea.       1.1 µm       100 µm PDMS (Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, ea.       6ray       1.1 µm       100 µm PDMS (Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Gray, a.       1.1 µm       100 µm PDMS (Polydimethylsiloxane)         PAL Smart SPME Arrow 110 mm: PA, Phase Thickness 100 um, Phase Length 20 mm, Gray, a.       1.1 µ	PAL Smart SPRE Arrow Method Development Kit       Light Rue       1.1 mm       120 µm Carbon-WR/PDMS (Carbon Wice Range/ Polydimethylsiloxane)       Highly volatile, 30–225 g/mol*         PAL Smart SPNE Arrow J.10 mm: Carbon-WR/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Light Blue, 3.p.k.       Light Blue       1.1 mm       120 µm Carbon-WR/PDMS (Carbon Wice Range/ Polydimethylsiloxane)       Highly volatile, 30–225 g/mol*         PAL Smart SPNE Arrow J.10 mm: DVB/Carbon-WR/PDMS, Phase Thickness 20 um, Phase Length 20 rm, Dark Gray, 9.       Dark Gray       1.1 mm       120 µm DVB/Carbon-WR/PDMS (Chron Wice Range/Polydimethylsiloxane)       Volatile and semivolatile, 40–275 g/mol*         PAL Smart SPNE Arrow J.10 mm: DVB/Carbon-WR/PDMS, Phase Thickness Length 20 mm, Volet, ea.       Volet       1.1 mm       120 µm DVB/Carbon-WR/PDMS (Divinylbenzene/ Polydimethylsiloxane)       Volatile and semivolatile, 40–275 g/mol*         PAL Smart SPNE Arrow J.10 mm: DVB/PDMS, Phase Thickness 120 um, Phase Length 20 mm, Volet, ea.       Volet       1.1 mm       120 µm DVB/PDMS (Divinylbenzene/ Polydimethylsiloxane)       Amines and polar compounds, 60-300 g/mol*         PAL Smart SPNE Arrow J.10 mm: PDMS, Phase Thickness 100 um, Phase Length 20 mm, Red, ea.       1.1 mm       100 µm PDMS (Polydimethylsiloxane)       Volatile, 60–275 g/mol* Polydimethylsiloxane)       Volatile, 60–275 g/mol* Polydimethylsilo

\*These molecular weight ranges are a reasonable approximation; however, end users should verify suitability for their specific application.



#### **Restek PAL SPME Arrow GC-Specific Conversion Kits**

Conversion kits designed for use with 1.1 and 1.5 mm Restek PAL SPME Arrows. Compatible with Smart SPME Arrows and non-smart SPME Arrows.

Due to the relatively large diameter of Restek PAL SPME Arrows, you must modify the GC inlet using an instrument-specific conversion kit prior to use. Converted inlets are compatible with all standard injection techniques (SPME Arrow, liquid syringe, headspace, etc.); so there is no need to switch inlets after installation.



Restek PAL SPME Arrow GC-Specific Conversion Kits are easy to install. For the weldment, install according to your instrument's owner's manual.

Description	Includes	Instrument	qty.	cat.#
Restek PAL SPME Arrow Conversion Kit	Topaz 1.8 mm ID straight/SPME inlet liner, 5-pk. (cat.# 23280); Ther- molite Plus septa, 3-pk. (cat.# 23864); Split/splitless weldment; large, canister-type filter (cat.# 27502); Septum nut for 6890 split/splitless weldments (cat.# 27503); Injector adaptor cup (cat.# 27496)	for Agilent 6890 Split/Splitless Injector (for canister-type filters)	kit	27492
Restek PAL SPME Arrow Conversion Kit with 1.1 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23280); 1.1 mm Microseal (cat.# 23232); Split/Splitless Weldment; Large Canister Type Filter (cat.# 27502); Adaptor Cup (cat.# 27496); Nut (cat.# 23228)	for Agilent 6890 Split/Splitless Injector (for canister-type filters)	kit	27356
Restek PAL SPME Arrow Conversion Kit vith 1.5 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23280); 1.5 mm Microseal (cat.# 23233); Split/Splitless Weldment and Large Canister Type Filter (cat.# 27502); Adaptor Cup (cat.# 27496); Nut (cat.# 23228)	for Agilent 6890 Split/Splitless Injector (for canister-type filters)	kit	27361
estek PAL SPME Arrow Conversion Kit	Topaz 1.8 mm ID straight/SPME inlet liner, 5-pk. (cat.# 23280); Thermo- lite Plus septa, 3-pk. (cat.# 23864); Agilent split/splitless weldment and septum nut (cat.# 27504); Injector adaptor cup (cat.# 27496)	for Agilent 7890 Split/Splitless Injector	kit	27493
estek PAL SPME Arrow Conversion Kit vith 1.1 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23280); 1.1 mm Microseal (cat.# 23232); Agilent Weldment (cat.# 27504); Adaptor Cup (cat.# 27496); Nut (cat.# 23228)	for Agilent 7890 Split/Splitless Injector	kit	27357
Restek PAL SPME Arrow Conversion Kit vith 1.5 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23280); 1.5 mm Microseal (cat.# 23233); Agilent Weldment (cat.# 27504); Adaptor Cup (cat.# 27496); Nut (cat.# 23228)	for Agilent 7890 Split/Splitless Injector	kit	27362
Restek PAL SPME Arrow Conversion Kit with 1.1 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23280); 1.1 mm Microseal (cat.# 23232); Agilent Weldment (cat.# 28933); Adaptor Cup (cat.# 27496); Nut (cat.# 23228)	for Agilent 8890 Split/Splitless Injector	kit	28934
testek PAL SPME Arrow Conversion Kit	Topaz 1.8 mm ID straight/SPME inlet liner, 5-pk. (cat.# 23279); Ther- molite Plus septa, 3 pk. (cat.# 23872); Injection port weldment (cat.# 27500); Needle guide/septum nut (cat.# 27501); Injector adaptor cup (cat.# 27497)	for Shimadzu GC-2010 Split/Splitless Injector (not compatible with SE or Plus models)	kit	27491
Restek PAL SPME Arrow Conversion Kit vith 1.1 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23279); 1.1 mm Microseal (cat.# 23232); Port Weldment (cat.# 27500 ); Adaptor Cup (cat.# 27497); Adaptor Kit (cat.# 23229)	for Shimadzu GC-2010 Split/Splitless Injector (not compatible with SE or Plus models)	kit	27355
Restek PAL SPME Arrow Conversion Kit vith 1.5 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23279); 1.5 mm Microseal (cat.# 23233); Port Weldment (cat.# 27500); Adaptor Cup (cat.# 27497); Adaptor Kit (cat.# 23229)	for Shimadzu GC-2010 Split/Splitless Injector (not compatible with SE or Plus models)	kit	27360
estek PAL SPME Arrow Conversion Kit	Topaz 1.8 mm ID straight/SPME inlet liner, 5-pk. (cat.# 23278); Premium nonstick BTO septa, 3-pk. (cat.# 27090); Septum cap (cat.# 27505); Lin- er cap/septum holder (cat.# 27506); Injector adaptor cup (cat.# 27498)	for Thermo TRACE 1300/1310, 1600/1610 Split/Splitless Injector	kit	27494
estek PAL SPME Arrow Conversion Kit vith 1.1 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23278); 1.1 mm Microseal (cat.# 23232); Liner Cap/Septum Holder (cat.# 27506); Adaptor Cup (cat.# 27498); Nut (cat.# 23230)	for Thermo TRACE 1300/1310, 1600/1610 Split/Splitless Injector	kit	27358
estek PAL SPME Arrow Conversion Kit vith 1.5 mm Merlin Microseal	Topaz 1.8 mm ID Straight/SPME inlet liner, 5-pk. (cat.# 23278); 1.5 mm Microseal (cat.# 23233); Liner Cap/Septum Holder (cat.# 27506); Adaptor Cup (cat.# 27498); Adaptor Kit (cat.# 23230)	for Thermo TRACE 1300/1310, 1600/1610 Split/Splitless Injector	kit	27363
estek PAL SPME Arrow Conversion Kit	2.0 mm ID straight inlet liner, 5-pk. (cat.# 22267); Premium nonstick BTO septa, 3-pk. (cat.# 27096); Septum holder and support (cat.# 27507); Liner cap (cat.# 27508); Injector adaptor cup (cat.# 27499)	for Thermo TRACE Ultra Split/Splitless Injector	kit	27495
estek PAL SPME Arrow Conversion Kit vith 1.1 mm Merlin Microseal	2.0 mm ID straight inlet liner, 5-pk. (cat.# 22267); 1.1 mm Microseal (cat.# 23232); Liner Cap (cat.# 27508); Adaptor Cup (cat.# 27499); Adaptor Kit (cat.# 23231)	for Thermo TRACE Ultra Split/Splitless Injector	kit	27359
Restek PAL SPME Arrow Conversion Kit vith 1.5 mm Merlin Microseal	2.0 mm ID straight inlet liner, 5-pk. (cat.# 22267); 1.5 mm Microseal (cat.# 23233); Liner Cap (cat.# 27508); Adaptor Cup (cat.# 27499); Adaptor Kit (cat.# 23231)	for Thermo TRACE Ultra Split/Splitless Injector	kit	27364




#### **Accessories for SPME Arrows**

Description	Instrument	qty.	cat.#
	for Agilent GC 6890/7890/8890 Split/Splitless Injector	ea.	27496
	for Shimadzu GC 2010 Split/Splitless Injector	ea.	27497
Injector adaptor cup	for Thermo GC TRACE 1300/1310, 1600/1610 Split/ Splitless Injector	ea.	27498
Injector adaptor cup Injection port weldment Needle guide/septum nut Split/splitless weldment; large, canis- ter-type filter Septum nut for split/splitless weldments Split/splitless weldment and septum nut Septum cap Liner cap/septum holder	for Thermo GC TRACE Ultra Split/Splitless Injector	ea.	27499
Injection port weldment	for Shimadzu GC 2010 Split/Splitless Injector	ea.	27500
Needle guide/septum nut	for Shimadzu GC 2010 Split/Splitless Injector	ea.	27501
Split/splitless weldment; large, canis- ter-type filter	for Agilent GC 6890 Split/Splitless Injector	ea.	27502
Septum nut for split/splitless weldments	for Agilent GC 6890/7890 Split/Splitless Injector	ea.	27503
Collis I collision and the contract of the con	for Agilent GC 7890 Split/Splitless Injector	ea.	27504
Split/splitless weldment and septum nut	for Agilent 8890 Split/Splitless Injector	ea.	28933
Septum cap	for Thermo GC TRACE 1300/1310, 1600/1610 Split/ Splitless Injector	ea.	27505
Liner cap/septum holder	for Thermo GC TRACE 1300/1310, 1600/1610 Split/ Splitless Injector	ea.	27506
Septum holder and support	for Thermo GC TRACE Ultra Split/Splitless Injector	ea.	27507
Liner cap	for Thermo GC TRACE Ultra Split/Splitless Injector	ea.	27508



#### More phase volume and robust construction mean Restek PAL SPME Arrows outperform traditional fibers.

- Rugged, stainless-steel construction ensures longer lifetimes.
- Faster extraction increases productivity.
- Better sensitivity for lower LODs.

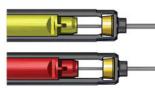



**1.5 mm Smart SPME Arrow** Surface area: 63 mm<sup>2</sup>, Volume: 12 μL

**1.1 mm Smart SPME Arrow** Surface: 44 mm<sup>2</sup>, Volume: 3.8 μL

**Smart SPME Fiber** Surface area: 9.4 mm², Volume: 0.6 μL



## **GC Inlet Liners for SPME**

RESTEK

RESTEK

#### Topaz 0.75 mm ID Straight/SPME Inlet Liner

for Bruker/Varian GCs w/1078/1079 Inlets

Geometry	Deactivation	Material	ID x OD x Length	qty	cat.#
Straight/SPME	Premium	Borosilicate Glass	0.75 mm x 5.0 mm x 54 mm	5-pk.	23465

#### Topaz 0.75 mm ID Straight/SPME Inlet Liner

for PerkinElmer Clarus 590/690 GCs

Geometry	Deactivation	Material	ID x OD x Length	qty	cat.#
Straight/SPME	Premium	Borosilicate Glass	0.75 mm x 6.35 mm x 78.5 mm	5-pk.	23803

#### Topaz 0.75 mm ID Straight/SPME Inlet Liner

for Agilent GCs

Geometry	Deactivation	Material	ID x OD x Length	qty	cat.#
Straight/SPME	Premium	Borosilicate Glass	0.75 mm x 6.35 mm x 78.5 mm	5-pk.	23434

#### Topaz 1.8 mm ID Straight/SPME Inlet Liner

for Shimadzu GCs 17A, 2010, 2014, and 2030 GCs

Geometry	Deactivation	Material	ID x OD x Length	qty	cat.#
Straight/SPME	Premium	Borosilicate Glass	1.8 mm x 5.0 mm x 95 mm	5-pk.	23279

#### Topaz 1.8 mm ID Straight/SPME Inlet Liner

for Agilent GCs

Geometry	Deactivation	Material	ID x OD x Length	qty	cat.#
Straight/SPME	Premium	Borosilicate Glass	1.8 mm x 6.5 mm x 78.5 mm	5-pk.	23280

#### Topaz 1.8 mm ID Straight/SPME Inlet Liner

for Thermo TRACE 1300/1310, 1600/1610 GCs w/SSL Inlets

Geometry	Deactivation	Material	ID x OD x Length	qty	cat.#
Straight/SPME	Premium	Borosilicate Glass	1.8 mm x 6.5 mm x 78.5 mm	5-pk.	23278

#### More GC inlet liners for SPME are available at www.restek.com

#### **SPME Performance Test Mix** (2 components)

- Essential mix for establishing the performance of SPME fibers and SPME Arrows.
- Verified composition and stability.

RESTEK

### Nitrobenzene (98-95-3)

2-Nitrotoluene (88-72-2)

Conc. in Solvent	CRM?	Min Shelf Life on Ship Date	Max Shelf Life on Ship Date	Shipping Conditions	Storage Temperature	qty.	cat.#
SPME Performance Test Mix							
1 μg/mL in water:methanol (99:1), 1 mL/ampul	Yes	6 months	36 months	Ambient	10 °C or colder	3-pk.	31015

BGB GC LC MS CE





RESTEK

RESTEK

## SPME Vials, Caps, and Septa

#### Magnetic Screw-Thread Caps, 18 mm

Description	Туре	Cap Size	Septa Material	qty.	cat.#
Magnetic Caps and Septa for	Screw-Thread	18-425	Blue PTFE/Silicone, 1.5 mm thick	100-pk.	23090
SPME	Screw-Thread	18-425	Blue PTFE/Silicone, 1.5 mm thick	1000-pk.	23091
	Screw-Thread	18-425	Red PTFE/Silicone, 1.9 mm thick	100-pk.	23092
Manual Constants	Screw-Thread	18-425	Red PTFE/Silicone, 1.9 mm thick	1000-pk.	23093
Magnetic Caps and Septa	Screw-Thread	18-425	PTFE/Red Chlorobutyl	100-pk.	23094
	Screw-Thread	18-425	PTFE/Red Chlorobutyl	1000-pk.	23095



23852

23854

#### SPME MicroCenter Caps and Septa

Description	Туре	Cap Size	Color	Septa Material	qty.	cat.#
	Screw-Thread	18-425		MicroCenter PTFE/Silicone, 0.040" (+/-0.005")	100-pk.	23852
	Screw-Thread	18-425		MicroCenter PTFE/Silicone, 0.040" (+/-0.005")	1000-pk.	23853
	Bi-Metal Crimp	20 mm	Blue	MicroCenter PTFE/Silicone, 0.065" (+/-0.005")	100-pk.	23854
CDME Viel Com	Bi-Metal Crimp	20 mm	Blue	MicroCenter PTFE/Silicone, 0.065" (+/-0.005")	1000-pk.	23855
SPME Vial Cap	Bi-Metal Crimp	20 mm	Red	MicroCenter PTFE/Silicone, 0.065" (+/-0.005")	100-pk.	23856
	Bi-Metal Crimp	20 mm	Red	MicroCenter PTFE/Silicone, 0.065" (+/-0.005")	1000-pk.	23857
	Steel Crimp	20 mm	Gold	MicroCenter PTFE/Silicone, 0.065" (+/-0.005")	100-pk.	23858
	Steel Crimp	20 mm	Gold	MicroCenter PTFE/Silicone, 0.065" (+/-0.005")	1000-pk.	23859
SPME Vial Septa,				MicroCenter PTFE/Silicone, 0.040" (+/-0.005")	100-pk.	23850
18 mm				MicroCenter PTFE/Silicone, 0.040" (+/-0.005")	1000-pk.	23851

Cat.# 23850 and 23851 not for use with 20 mm caps.

#### Headspace Crimp Vials, 20 mm

Modification	Туре	Volume	Color	Deactivation	Size	qty.	cat.#
Flat Bottom	20 mm Crimp-Top	6 mL	Clear		22 x 38 mm	100-pk.	21166
Flat Bottom	20 mm Crimp-Top	6 mL	Clear		22 x 38 mm	1000-pk.	21167
Flat Bottom	20 mm Crimp-Top	10 mL	Clear		23 x 46 mm	100-pk.	24683
Flat Bottom	20 mm Crimp-Top	10 mL	Clear		23 x 46 mm	1000-pk.	24684
Rounded Bottom	20 mm Crimp-Top	10 mL	Clear		23 x 46 mm	100-pk.	21164
Rounded Bottom	20 mm Crimp-Top	10 mL	Clear		23 x 46 mm	1000-pk.	21165
Rounded Bottom	20 mm Crimp-Top	10 mL	Clear	Deactivated	23 x 46 mm	1000-pk.	21165-221
Flat Bottom	20 mm Crimp-Top	20 mL	Clear		23 x 75 mm	100-pk.	24685
Flat Bottom	20 mm Crimp-Top	20 mL	Clear		23 x 75 mm	1000-pk.	24686
Rounded Bottom	20 mm Crimp-Top	20 mL	Clear		23 x 75 mm	100-pk.	21162
Rounded Bottom	20 mm Crimp-Top	20 mL	Clear		23 x 75 mm	1000-pk.	21163
Flat Bottom	20 mm Crimp-Top	27 mL	Clear		30 x 60 mm	100-pk.	21160
Flat Bottom	20 mm Crimp-Top	27 mL	Clear		30 x 60 mm	1000-pk.	21161
	Flat Bottom Flat Bottom Flat Bottom Rounded Bottom Rounded Bottom Flat Bottom Flat Bottom Rounded Bottom Rounded Bottom Flat Bottom	Flat Bottom     20 mm Crimp-Top       Rounded Bottom     20 mm Crimp-Top       Flat Bottom     20 mm Crimp-Top       Flat Bottom     20 mm Crimp-Top       Flat Bottom     20 mm Crimp-Top       Rounded Bottom     20 mm Crimp-Top       Rounded Bottom     20 mm Crimp-Top       Flat Bottom     20 mm Crimp-Top       Rounded Bottom     20 mm Crimp-Top       Rounded Bottom     20 mm Crimp-Top       Flat Bottom     20 mm Crimp-Top       Flat Bottom     20 mm Crimp-Top	Flat Bottom     20 mm Crimp-Top     6 mL       Flat Bottom     20 nm Crimp-Top     6 mL       Flat Bottom     20 nm Crimp-Top     10 mL       Flat Bottom     20 nm Crimp-Top     10 mL       Flat Bottom     20 nm Crimp-Top     10 mL       Rounded Bottom     20 nm Crimp-Top     10 mL       Rounded Bottom     20 nm Crimp-Top     10 mL       Rounded Bottom     20 nm Crimp-Top     10 mL       Flat Bottom     20 nm Crimp-Top     20 mL       Flat Bottom     20 nm Crimp-Top     20 mL       Flat Bottom     20 nm Crimp-Top     20 mL       Rounded Bottom     20 nm Crimp-Top     20 mL       Flat Bottom     20 nm Crimp-Top     20 mL       Flat Bottom     20 nm Crimp-Top     20 mL       Rounded Bottom     20 nm Crimp-Top     20 mL       Flat Bottom     20 nm Crimp-Top     20 mL       Rounded Bottom     20 nm Crimp-Top     20 mL	Flat Bottom20 mm Crimp-Top6 mLClearFlat Bottom20 mm Crimp-Top6 mLClearFlat Bottom20 mm Crimp-Top10 mLClearFlat Bottom20 mm Crimp-Top10 mLClearFlat Bottom20 mm Crimp-Top10 mLClearRounded Bottom20 mm Crimp-Top10 mLClearRounded Bottom20 mm Crimp-Top10 mLClearRounded Bottom20 mm Crimp-Top10 mLClearFlat Bottom20 mm Crimp-Top20 mLClearFlat Bottom20 mm Crimp-Top20 mLClearRounded Bottom20 mm Crimp-Top20 mLClearFlat Bottom20 mm Crimp-Top27 mLClear	Flat Bottom       20 mm Crimp-Top       6 mL       Clear         Flat Bottom       20 mm Crimp-Top       6 mL       Clear         Flat Bottom       20 mm Crimp-Top       10 mL       Clear         Flat Bottom       20 mm Crimp-Top       10 mL       Clear         Flat Bottom       20 mm Crimp-Top       10 mL       Clear         Rounded Bottom       20 mm Crimp-Top       10 mL       Clear         Rounded Bottom       20 mm Crimp-Top       10 mL       Clear         Rounded Bottom       20 mm Crimp-Top       10 mL       Clear         Flat Bottom       20 mm Crimp-Top       10 mL       Clear         Flat Bottom       20 mm Crimp-Top       20 mL       Clear         Flat Bottom       20 mm Crimp-Top       20 mL       Clear         Rounded Bottom       20 mm Crimp-Top       20 mL       Clear         Flat Bottom       20 mm Crimp-Top       20 mL       Clear	Flat Bottom         20 mm Crimp-Top         6 mL         Clear         22 x 38 mm           Flat Bottom         20 mm Crimp-Top         6 mL         Clear         22 x 38 mm           Flat Bottom         20 mm Crimp-Top         6 mL         Clear         22 x 38 mm           Flat Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm           Flat Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm           Rounded Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm           Rounded Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm           Rounded Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm           Rounded Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm           Flat Bottom         20 mm Crimp-Top         20 mL         Clear         23 x 46 mm           Flat Bottom         20 mm Crimp-Top         20 mL         Clear         23 x 45 mm           Flat Bottom         20 mm Crimp-Top         20 mL         Clear         23 x 75 mm           Rounded Bottom         20 mm Crimp-Top         20 mL         Clear         23 x 75 mm	Flat Bottom         20 mm Crimp-Top         6 mL         Clear         22 x 38 mm         100-pk.           Flat Bottom         20 mm Crimp-Top         6 mL         Clear         22 x 38 mm         100-pk.           Flat Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm         100-pk.           Flat Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm         100-pk.           Flat Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm         100-pk.           Rounded Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm         100-pk.           Rounded Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm         100-pk.           Rounded Bottom         20 mm Crimp-Top         10 mL         Clear         23 x 46 mm         100-pk.           Rounded Bottom         20 mm Crimp-Top         20 mL         Clear         23 x 46 mm         100-pk.           Flat Bottom         20 mm Crimp-Top         20 mL         Clear         23 x 75 mm         100-pk.           Rounded Bottom         20 mm Crimp-Top         20 mL         Clear         23 x 75 mm         100-pk.           Rounded Bottom <t< td=""></t<>

Vial-to-instrument compatibility is designated in Instrument Reference Chart for Headspace Vials at www.restek.com.

#### Headspace Screw-Thread Vials, 18 mm

Description	Modification	Туре	Volume	Color	Size	qty.	cat.#
	Rounded Bottom	18-425 Screw-Thread	20 mL	Clear	22 x 75 mm	100-pk.	23082
	Rounded Bottom	18-425 Screw-Thread	20 mL	Clear	22 x 75 mm	1000-pk.	23083
	Rounded Bottom	18-425 Screw-Thread	20 mL	Amber	22 x 75 mm	100-pk.	23086
Headspace Vial	Rounded Bottom	18-425 Screw-Thread	20 mL	Amber	22 x 75 mm	1000-pk.	23087
Rounded Bottom	Rounded Bottom	18-425 Screw-Thread	10 mL	Clear	22 x 45 mm	100-pk.	23084
	Rounded Bottom	18-425 Screw-Thread	10 mL	Clear	22 x 45 mm	1000-pk.	23085
	Rounded Bottom	18-425 Screw-Thread	10 mL	Amber	22 x 45 mm	100-pk.	23088
	Rounded Bottom	18-425 Screw-Thread	10 mL	Amber	22 x 45 mm	1000-pk.	23089

Caps not included (sold separately).



21166



23082





## **Merlin Microseals for SPME Arrow**

#### Merlin Microseal Nut for SPME Arrow for Agilent GCs

Description	Includes	Instrument	Туре	Vendor cat.#	qty.	cat.#
Merlin Microseal Nut for SPME Arrow	nut (1)	Agilent	SPME Arrow Application Nut (3 to 100 psi)	1000AG	ea.	23228

#### Merlin Microseal Adaptor Kit for SPME Arrow for Shimadzu 2010, 2025, and 2030 GCs

Description	Includes	Instrument	Туре	Vendor cat.#	qty.	cat.#
Merlin Microseal Adaptor Kit for SPME Arrow	nut (1); adaptor (1); O-rings (4)	Shimadzu 2010, 2025, and 2030	SPME Arrow Application Adaptor Kit (3 to 100 psi)	1000SH	kit	23229

#### Merlin Microseal Nut for SPME Arrow for Thermo TRACE 1300/1310, 1600/1610 GCs

Description	Includes	Instrument	Туре	Vendor cat.#	qty.	cat.#
Merlin Microseal Nut for SPME Arrow	nut (1)	TRACE 1300/1310  TRACE 1600/1610	SPME Arrow Application Nut (3 to 100 psi)	1000TS	ea.	23230

#### Merlin Microseal Adaptor Kit for SPME Arrow for Thermo TRACE Ultra GCs

Description	Includes	Instrument	Туре	Vendor cat.#	qty.	cat.#
Merlin Microseal Adaptor Kit for SPME Arrow	nut (1); adaptors (3); O-rings (4)	Thermo TRACE Ultra	SPME Arrow Application Adaptor Kit (3 to 100 psi)	1000TU	kit	23231



#### Description Type

**Merlin Microseal Replacement Septa** 

		cat.m		
	General-Purpose Microseal (most applications, 3 to 100 psi)	410	ea.	22812
	Low-Pressure Microseal (1 to 45 psi)	310	ea.	22815
Deale com out Microsof	Microseal for Traditional SPME Fiber Applications (3 to 100 psi)	21-01	ea.	22782
Replacement Microseal	Microseal for 1.1 mm SPME Arrow Applications (3 to 100 psi)	1100	ea.	23232
-	Microseal for 1.5 mm SPME Arrow Applications (3 to 100 psi)	1500	ea.	23233
	Microseal for 26 gauge or 23/26 gauge tapered needles (5 to 100 psi)	610	ea.	22264

## Our product line is continually expanding! See what's new at www.restek.com/SPME

Vendor

cat #

qty.

cat.#

BGB	sclic	
DGD	MS CE DR-BGB-2404	www.bgb-shop.com
Switzerland Austria	BGB Analytik AG • Rohrmattstrasse 4 • 4461 Böckten • Phone +41 61 991 00 46 • Fax +41 61 991 00 25 • sales⊚ BGB Analytik Vertrieb GmbH • Mühlestraße 1 • 79539 Lörrach • Phone +49 7621 5884270 • Fax +49 7621 58842	, , , , , , , , , , , , , , , , , , ,
Turkey and Azerbaijan	BGB Anlaytik Vertreb Ginbi i * Muniestralse i * 79339 Lonach * Filone +49 7021 3664270 * Fax +49 7021 36642 BGB Analytik Laboratuvar Malzemeleri Ticaret LTD.ŞTİ • Starport Residence • Yenişehir Mah. Osmanlı Bulvarı. • 3 34912 Kurtköy, Pendik, İstanbul • Phone +90 2169092048 • sales.tr@bgb-analytik.com	<b>3 3 7</b>

